
Q-smart HE

Compact High-Energy pulsed Nd:YAG lasers with excellent beam quality and versatility



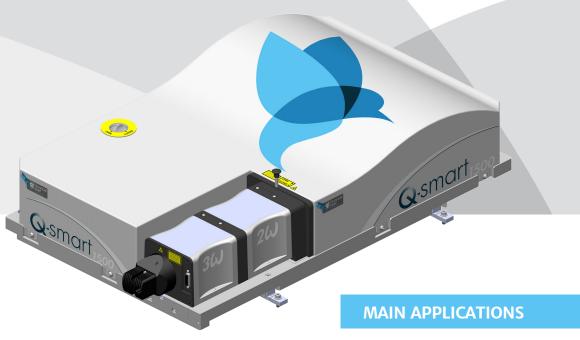


Up to 2.3 J @ 1064 nm

MAIN FEATURES

- Robust and field proven technology
- Built to last thanks to ceramic reflectors and long flashlamp life time warranty
- Plug & play harmonic modules with automatic phase-matching
- Cables and cooling lines fully disconnectable
- · Easy to use and maintain
- Optional chiller
- · Intuitive GUI interface
- SLM option (Single Longitudinal Mode)

- LiDAR
- INSTRUMENTATION
- PLD
- · DYE, OPO & Ti:Sa PUMPING
- SPECTROSCOPY
- · LIF
- COMBUSTION


Typical beam profiles

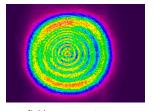

Near field 1.5 J @ 1064 nm, 10 Hz



Far field 1.5 J @ 1064 nm, 10 Hz

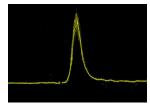

6 ns typical temporal profile @ 1064 nm (1 GHz oscilloscope)

www.quantel-laser.com

Please contact Lumibird to find the best match for your needs and compatibility between options.





Q-smart HE

Compact High-Energy pulsed Nd:YAG lasers with excellent beam quality and versatility



SPECIFICATIONS

		Q-smart 1200		Q-smart 1500		Q-smart 2300
Repetition rate (Hz)		10	20	10	20	10
Energy per pulse (mJ)	1064 nm	1200	1000	1500	1400	2300
	532 nm	575 / 650 ⁽¹⁾	480 / 540 (1)	750 / 850 (1)	660 / 750 ⁽¹⁾	1150/1300 ⁽¹⁾
	355 nm	280 / 350 (2)	220 / 270 (2)	400 / 520 (2)	360 / 460 (2)	620 / 850 ⁽²⁾
	266 nm	110	80	130	100	200
Pulse duration (ns) (3)	1064 nm	5 - 10				
Beam diameter (mm) (4)	1064 nm	≤ 10 ≤			≤ 12	
Beam divergence (mrad) (5)	1064 nm	≤ 0.5				
M ^{2 (6)}	1064 nm	≤2				
Spatial profile @ 1064 nm ⁽⁷⁾ (fit to Gaussian)	Near field (8)	≥ 0.7				
	Far field (9)	≥ 0.9				
Polarization ratio (%) (10)	1064 nm	≥ 90 ≥ 80				

- (1) 532 nm high energy version
- (2) 355 nm high energy version
- (3) Measured at FWHM with fast photodiode and 1 GHz oscilloscope
- (4) At the output of the laser
- (5) Full angle, at 1/e² of the peak
- (6) At 1/e2 of the peak, measured by Spricon LBA FWB
- (7) Least square fit to Gaussian (perfect fit = 1)
- (8) Measured at 1 m from laser output
- (9) Measured at focal plane of a 2 m focus lens
- (10) Polarization is horizontal @ 1064, 355 & 266 nm and vertical @ 532 nm


1064 nm	± 2 (0.6)
532 nm	± 4 (1.3)
355 nm	± 6 (2)
266 nm	± 8 (2.6)
1064 nm	± 3
532 nm	± 5
355 nm	± 5
266 nm	± 10
1064 nm	< 40
Standard (14)	≤ 0.7
SLM (15) option	≤ 0.005
	532 nm 355 nm 266 nm 1064 nm 532 nm 355 nm 266 nm 1064 nm

- (11) Peak-to-peak (RMS), 100% of shots
- (12) Over 8 hours for $\Delta T^{\circ} \leq \pm 3^{\circ}C$
- (13) Measured with Spiricon LBA-100, rms, on 200 pulses at the focal plane
- of a 2 m focus lens
- (14) Measured at FWHM with a grating spectrometer with 0.045 cm⁻¹ resolution
- (15) Measured at FWHM with a slow scan Fabry-Perot etalon, ≤ 20% energy reduction @ 1064 nm



OTHER INFORMATION					
Power requir	ements	240 V (110 V on request) 16 to 20 A, 50/60 Hz			
Cooling		Water to water Optional chiller (stand alone or 19» rack)			
Operating ter	mperature	+ 18 °C to + 28 °C			
Storage temp	perature (16)	- 10 °C to + 50 °C			
Cable length	(m)	3 (17)			
Flashlamps v	varranty	50 million shots (18)			
Weight (kg)	Laser head	45			
	Harmonic modules	2.1			
	Integrated cooling & electronics	50 / 70 (19)			

- (16) System rinsed and drained with ethylene glycol/water mixture (17) Other lengths on request. Some losses are to be expected.
- (18) 80% of energy, or 1 year, whichever comes first (19) Q-smart 1200-1500 / Q-smart 2300

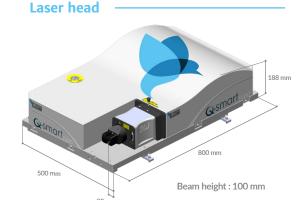

Integrated cooling & electronics



www.quantel-laser.com



